

# Delta license key申请步骤

1. 打开网址：<https://nmrsupport.jeol.com>
2. 申请登录账户：

JEOL Delta / NMR Support Site

LOGIN | REGISTER

JEOLUSA.com | JEOL Resonance

JEOL Solutions for Innovation

Input your search...

Home

## Welcome to JEOL Delta / NMR Support

The JEOL Delta / NMR support web site provides support for users interested in the Delta NMR Software package and JEOL's NMR Spectrometers. The web site includes Delta NMR software kits for NMR data processing only and kits for JEOL NMR Spectrometer control with data processing. Areas for Delta License Key generation, discussion forums, FAQ's, and documentation downloads are also available on this web site.

**To get started, simply:**

- 1. Open a User Account**  
Registration to this site is moderated - to gain full access, be sure to **also submit your spectrometer information** so that we can assign you the proper authorization for software downloads and/or license key generation.
- 2. Login to Your Existing Account**  
You will then see content and features available for your account type.
- 3. Migrate Your Account from AI-SUP, the old Delta support web site**  
If you are a user migrating from [ai-sup.jeol.com](https://ai-sup.jeol.com) please **click to login and update your account and spectrometer information in order to gain access to the appropriate area(s) of the site.**

Once your account registration and spectrometer information is reviewed, you will be granted the appropriate access to other areas of the site.

### 3. 填写账户信息：（\*内容为必填项）

JEOL Delta / NMR Support Site LOGIN | REGISTER

JEOLUSA.com | JEOL Resonance f t in YouTube

**JEOL**  
Solutions for Innovation

Input your search...

Home

## JEOL Delta/NMR Support Site Account Registration, Page 1

Completely fill out this form. Your request will be forwarded to the site administrator. You will be contacted once your account has been approved.

**Prefix:**

**\* First {Given} Name:**

**姓名**

**\* Family {Last} Name:**

**名**

**Suffix:**

**Display Name:**

**职业身份**

**\* Job Title:**

**公司/学校名称** **\* Company, or University Name, No Abbreviations Please:**

**Department:**

**邮箱地址** **\* Email Address:**

**邮箱再次确认** **\* Email Address (confirm):**

How did you find out about Delta?

**Make a Selection...**

- Faculty
- Post Doctoral Fellow
- Graduate Student
- Student
- NMR Technical Staff**
- Research Scientist
- Research Staff
- IT Staff
- Other
- JEOL Technical Staff
- JEOL Service Staff
- JEOL Sales Staff

### 3. 填写账户信息:

	<b>* Email Address:</b> <input type="text"/>
	<b>* Email Address (confirm):</b> <input type="text"/>
	<b>How did you find out about Delta:</b> JEOL NMR Staff ▼
接受使用条款	<b>* Acceptance of Terms of Use:</b> <input checked="" type="checkbox"/> <a href="#">Terms of Use</a>
地址信息	<b>* Address:</b> <input type="text" value="East Zhongguancun Road"/>
	<b>Unit:</b> <input type="text"/>
	<b>Address (2):</b> <input type="text"/>
	<b>Address (3):</b> <input type="text"/>
	<b>* City:</b> <input type="text" value="Beijing"/>
	<b>* Country:</b> <input type="text" value="China"/>
	<b>* State / Region:</b> <input type="text" value="Beijing"/> ▼
	<b>* Postal Code:</b> <input type="text" value="100224"/>
电话	<b>* Telephone:</b> <input type="text"/>
	<b>Cellular Phone:</b> <input type="text"/>
	<b>Fax:</b> <input type="text"/>
	<b>Time Zone:</b> (UTC-12:00) International D ▼
用户名	<b>* Username:</b> <input type="text"/>
密码	<b>Password:</b> <input type="text"/>
确认密码	<b>Confirm Password:</b> <input type="text"/>

#### 4. 点 “Continue”继续完善后面的仪器信息

Spectrometer Brand	JEOL
JEOL Spectrometer Type	ECZR
Spectrometer Field	600
Workstation	Intel PC
Operating System	Windows
Operating System Version	7 32 bit

Do you need Delta for:  Data Processing Only  
 Data Processing and JEOL NMR Machine Control (JEOL NMR Spectrometers Only)

JEOL Spectrometer serial number (format 'NM1xxxxxx' on back of main unit)  
Multiple JEOL NMR Spectrometer Laboratories only need to register one JEOL NMR system

Are you a JEOL employee or representative?  No  Yes


Submit

均需勾选

仪器序列号，只是申请数据处理的可以不填写

5. 填写完成点“submit”提交

6. 等待邮件通知，注册完成之后，返回网页(<https://nmrsupport.jeol.com>)登录

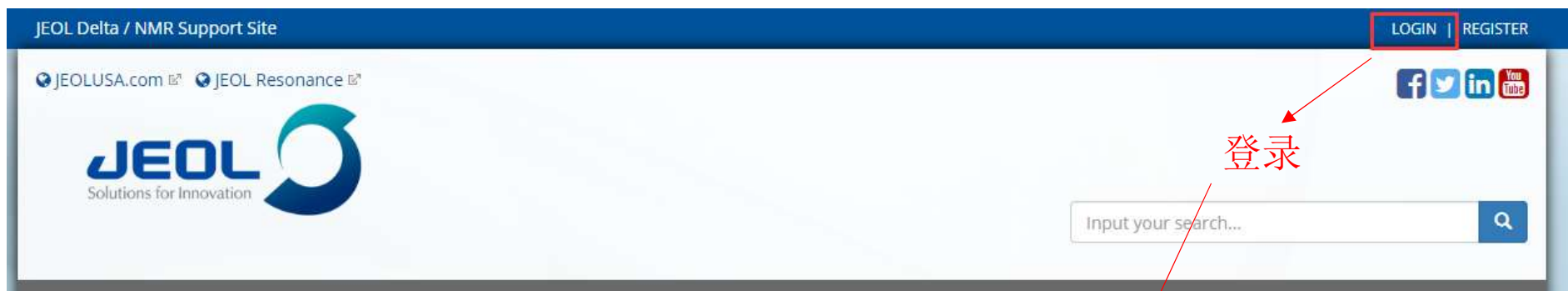

## 7. 填写计算机信息

CAREFULLY READ THE FOLLOWING TERMS AND CONDITIONS. YOU AGREE TO ALL TERMS AND CONDITIONS BY INSTALLING THE DELTA NMR SOFTWARE PACKAGE OR USING ANY PART THERE OF.

License: With the payment of any and all applicable License Fees upon the delivery of Delta NMR Software ('Software'), JEOL USA, INC. ('Licensor') grants to the Licensee a License to use without exclusivity a single non-transferable, non-sublicensable copy of the Software on a single CPU. No title to, ownership of, or rights to the software is hereby transferred to the licensee. The Software and its ancillary documentation may not be modified, reverse engineered, sublicensed, distributed, sold as a service on a time-sharing system, or used for any use other than the licensee's own internal use. This Software is composed of media, computer programs, and supporting documentation. In all cases the Software and its associated documentation may not be copied in whole or in part by any means or in any way disclosed to a third party. Upon License termination all Media, Software and Documentation must be removed from the CPU's and premises of the Licensee and returned to the licensor. The obligations of non-replication and non-disclosure set forth above shall survive any termination of this license.

Accept Terms  Reject Terms

接受使用条款

PackageType  Data Processing Only

Full Key with Instrument Control

根据邮件中的审核结果选择安装类型

Workstation Intel PC


Operating System Windows

Operating System Version 7 32 bit

Delta Machine ID UN - NI - NC - AT - HW

点击软件中的Options\_License Key获得软件ID

NOTE: The Delta Machine ID is generated by Delta not the operating system. It is not the MAC address.

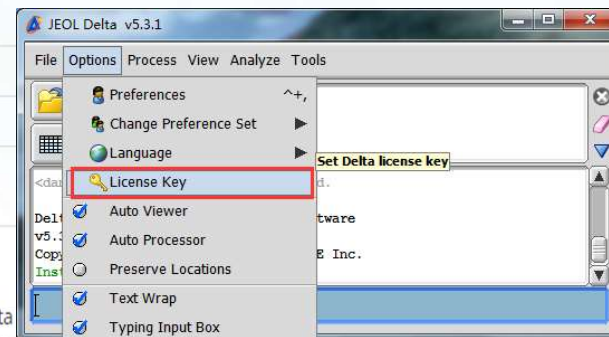

Submit

提交，获得license key

